


OM-MOUNT PHOTOMICRO ADAPTER L

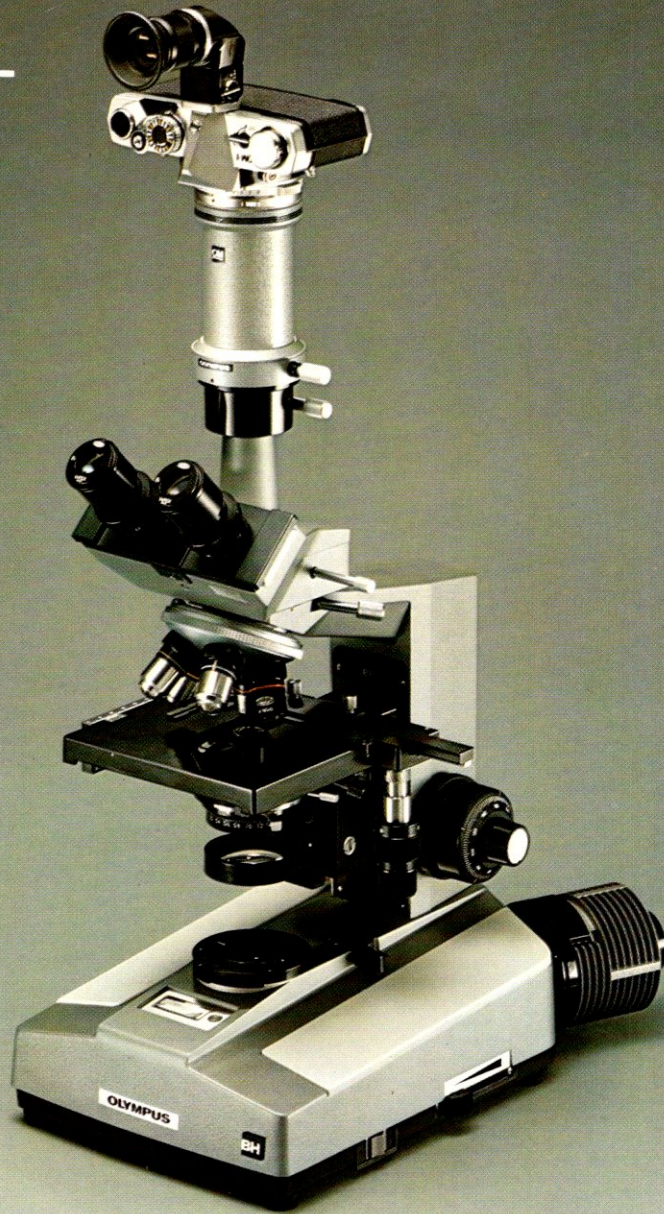

Summary

- A simple photomicrographic adapter for use of OM camera bodies in conjunction with microscopes.

Main Characteristics

- A special eyepiece adapter is required to mount an eyepiece into the photo tube of an ordinary microscope for objective magnifications up to 20X.
- No special eyepiece adapter is required to mount an eyepiece to the Olympus Research Microscope VANOX or Photomicrographic Stand (p.612). It assures sharply focused photomicrographs over the whole range of magnifications, eliminating any shutter vibration.

Notes • Related Units

- A special eyepiece adapter (p.611) is required to mount an eyepiece to an ordinary microscope.
- If used with the Olympus Research Microscope VANOX or Photomicrographic Stand PM-PSS (p.612), no special eyepiece adapter is required.
- Use the focusing screen 1-12 for photomicrography (p.201).
- The Varimagni Finder (p.212) is recommended in conjunction with the Eyecoupler (p.221) for sharp focusing in photomicrography.

OLYMPUS

OM SYSTEM OM-MOUNT PHOTOMICRO ADAPTER L

● Meter Coupling

Sets the exposure meter built into the camera body to the exposure Value necessary for photo micrography.

● Camera Mounting Ring

Connects to the OM body by means of bayonet mounting in the same manner as normal lenses.

● Microscope Mounting Ring

Fit the Adapter L on the eyepiece adapter, Research Microscope VANOX or Photomicrographic Stand PM-PSS.




● Adapter Release Button

Press the button and rotate the adapter L counter-clockwise by 70° to remove. For mounting, do not press the button.

● Microscope Clamping Screw

Clamp the adapter with screw provided.

Actual size

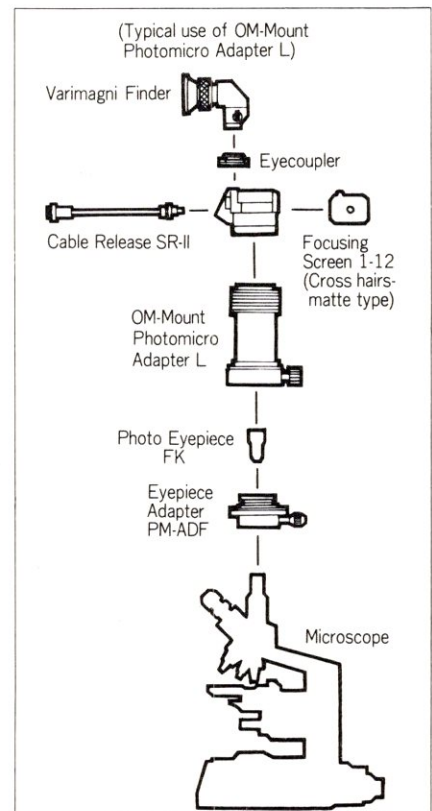

OLYMPUS